



Improved purification of histidine-tagged proteins with Ni Sepharose high performance

Intellectual Property Notice: The Biopharma business of GE Healthcare was acquired by Danaher on 31 March 2020 and now operates under the Cytiva™ brand. Certain collateral materials (such as application notes, scientific posters, and white papers) were created prior to the Danaher acquisition and contain various GE owned trademarks and font designs. In order to maintain the familiarity of those materials for long-serving customers and to preserve the integrity of those scientific documents, those GE owned trademarks and font designs remain in place, it being specifically acknowledged by Danaher and the Cytiva business that GE owns such GE trademarks and font designs.

cytiva.com

GE and the GE Monogram are trademarks of General Electric Company.
Other trademarks listed as being owned by General Electric Company contained in materials that pre-date the Danaher acquisition and relate to products within Cytiva's portfolio are now trademarks of Global Life Sciences Solutions USA LLC or an affiliate doing business as Cytiva.
Cytiva and the Drop logo are trademarks of Global Life Sciences IP Holdco LLC or an affiliate.
All other third-party trademarks are the property of their respective owners.
© 2020 Cytiva
All goods and services are sold subject to the terms and conditions of sale of the supplying company operating within the Cytiva business. A copy of those terms and conditions is available on request. Contact your local Cytiva representative for the most current information.
For local office contact information, visit [cytiva.com/contact](https://www.cytiva.com/contact)

Improved Purification of Histidine-tagged Proteins with Ni Sepharose High Performance

Lars C. Andersson[†], Lars Anderson*, Ann Bergh[†], Anna Heijbel[†], Helena Lindgren[†], Jon Lundqvist[†], Henrik Stålbbrand*, Karin Torstenson[†] and Katarina Öberg[†].

[†] GE Healthcare Bio-Sciences AB, Björkgatan 30, SE-751 84 Uppsala, Sweden

* Dept. Biochemistry, Lund University, P.O. Box 124, SE-221 00 Lund, Sweden

Introduction

The (histidine)₆-tag has become the most used affinity tag due to its small size, strong metal ion binding, and ability to bind under denaturing conditions. Immobilized Metal Ion Affinity Chromatography (IMAC) with Ni²⁺ is the standard method for purifying histidine-tagged proteins. Ni Sepharose™ High Performance medium is optimized for purification of such proteins.

Ni Sepharose High Performance

- Is precharged with Ni²⁺ and available in different formats, including prepacked 1 ml and 5 ml HisTrap™ HP columns.
- Consists of 34 µm beads. The small bead-size gives narrow peaks with high target protein concentration.
- Is compatible with a wide range of buffers and additives, including denaturants, detergents, and reducing agents.
- Has very low nickel ion leakage.
- Has very high protein binding capacity.

For detailed information see Data File 18-1174-40.

Conclusions

The following results were obtained with Ni Sepharose High Performance:

- Low leakage of Ni²⁺ was confirmed. A low leakage minimizes potential problems such as Ni²⁺-induced oligomerization and precipitation of target protein, and loss of binding capacity.
- The excellent binding properties of the medium allowed high flowrate purifications.
- A number of histidine-tagged proteins from different sources, with different M_r and with different expression levels were easily purified.
- 60 mg of purified protein could be obtained in a single run on a 1 ml column.



Material and Methods

Chromatography conditions were, unless otherwise stated:

Medium:	Ni Sepharose High Performance prepacked in HisTrap HP 1 ml columns or packed in 5x50 mm glass columns.
Samples: Clarified	Cell extracts from <i>E. coli</i> or <i>Pichia. pastoris</i> . by centrifugation and filtration; containing 0.5 M NaCl and imidazole at a concentration appropriate for each target protein.
Flow rate:	1 ml/min.
Binding buffer:	20 mM sodium phosphate, 0.5 M NaCl, x mM imidazole, pH 7.4.
Optimization:	The optimal imidazole concentration, to obtain the best purity and yield, differs from protein to protein. A small number of screening runs, including a run with an imidazole gradient from 5 mM, will facilitate finding a suitable imidazole concentration for optimal results.
Elution buffer:	20 mM sodium phosphate, 0.5 M NaCl, 500 mM imidazole, pH 7.4.

Chromatography was performed on ÄKTAexplorer™ 10 or 100 systems.

Elution was performed either with a step or a linear imidazole gradient.

Fractions of 1 ml were collected. SDS-electrophoresis was performed with ExcelGel™ SDS Gradient 8–18 gels.


Low nickel ion leakage

A. Comparison of Ni²⁺ leakage from three IMAC media at pH 4

- 10 column volumes of pH 4.0 buffer (no protein).
- Ni²⁺ still bound was eluted with EDTA and the amount was determined spectrophotometrically (372 nm).

Results

- Ni²⁺ leakage was considerably lower for Ni Sepharose High Performance than for the other media.

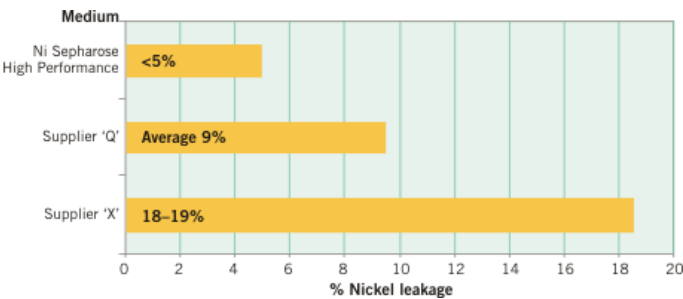

B. Ni²⁺ concentration in purified target protein pools from two different IMAC media

- Samples: Histidine-tagged proteins in *E. coli* and *P. pastoris* cell extracts.
- IMAC: Equal volumes of identical extracts applied to two different IMAC media. 5 mM imidazole for equilibration, sample application, and wash. Elution with linear imidazole gradient (chromatograms not shown). Pool volumes differed due to different peak widths.
- Ni²⁺ concentration was determined by atomic absorption spectrophotometry in eluted pools of target protein (of very similar purity).

Results

- Lower Ni²⁺ content in pools from Ni Sepharose High Performance.

Note: Potential problems due to Ni²⁺ contamination may vary with the molar ratio Ni²⁺/protein. The low ratio obtained for Ni Sepharose High Performance is due to the low Ni²⁺ concentration and the high target protein concentration obtained with small (34 µm) beads.

Histidine-tagged green fluorescent protein from *E. coli*. Data from final pool of purified protein:

	Ni Sepharose HP	Supplier 'Q'
volume of final pool	7 ml	11 ml
µg Ni ²⁺ /L (=ppb)	185	231
total amount of Ni ²⁺ in pool (nmol)	22	43
protein conc (mg/ml)	1.60	0.91
molar ratio Ni ²⁺ /protein	0.055	0.121



Histidine-tagged hydrolase from *P. pastoris*. Data from final pool of purified protein:

	Ni Sepharose HP	Supplier 'Q'
volume of final pool	5 ml	8 ml
µg Ni ²⁺ /L (=ppb)	87	129
total amount of Ni ²⁺ in pool (nmol)	7	18
protein conc (mg/ml)	n.d.	n.d.
molar ratio Ni ²⁺ /protein	n.d	n.d

n.d. = not determined

Purifications at high flow rates

1, 2 or 4 ml/min on 1 ml HisTrap HP columns

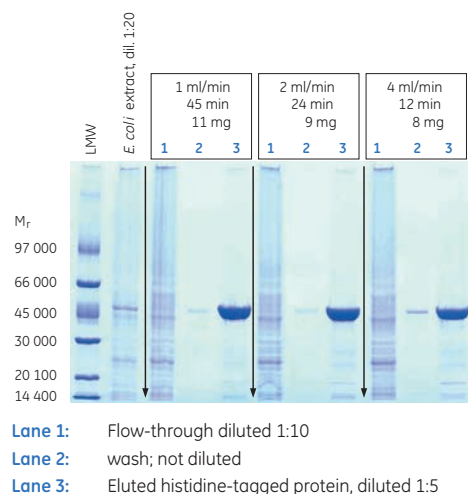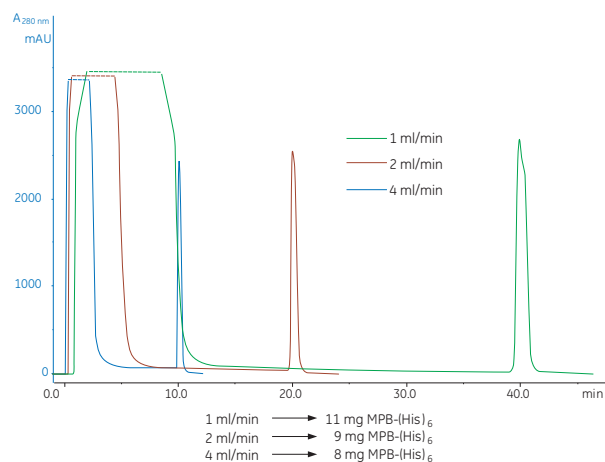

Results

- Samples: *E. coli* extract with histidine-tagged maltose binding protein (MBP-(His)₆, $M_r \sim 44\ 000$) at ~ 1.7 mg/ml.
- IMAC: At 1, 2 or 4 ml/min on 1 ml HisTrap HP. 8 ml extract applied (13.6 mg MBP-(His)₆). 35 mM imidazole during equilibration, sample application, and wash (see Optimization in Materials and Methods). Step elution with 500 mM imidazole.
- Only a slight decrease in yield was seen in the high-flow runs.
- SDS-PAGE showed very similar purity at low and high flow rates.

Purification of a high-MW histidine-tagged mannanase expressed in *E. coli*

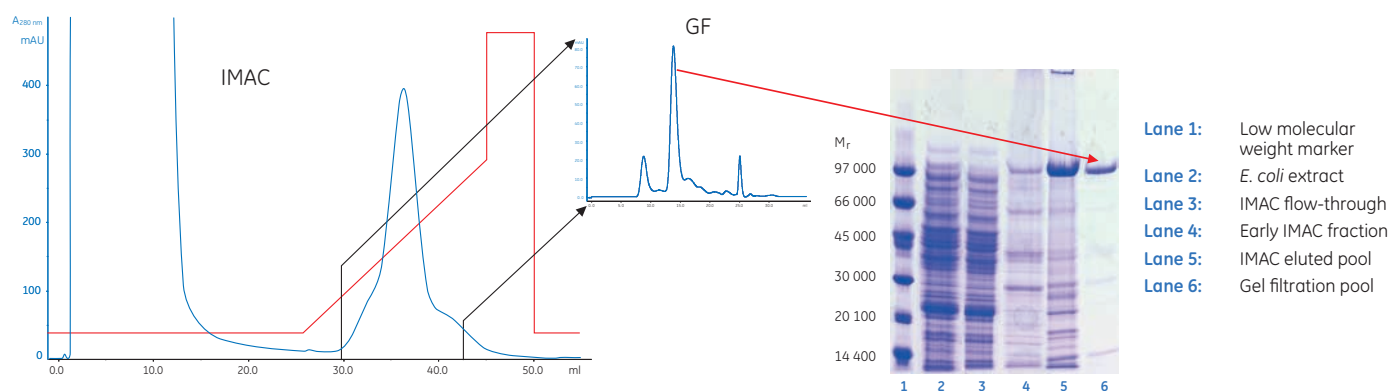


Fig 4. Purification by IMAC and gel filtration. Black lines in IMAC chromatogram indicate the IMAC pool taken to gel filtration.

Results

- Sample: *E. coli* extract with low-level expression of a histidine-tagged mannanase, Man 26A, from *Cellulomonas fimi* ($M_r \sim 100\ 000$).
 - IMAC: 10 ml extract applied to a 1 ml column with Ni Sepharose High Performance. 30 mM imidazole during equilibration, sample application, and wash (see Optimization in Materials and Methods). Elution with a 25 ml linear imidazole gradient to 300 mM.
 - Gel filtration: Superdex™ 200 10/300 GL.
 - Enzymatically active purified Man 26A.
- Note:* Several contaminants were co-purified with the full-length target protein. Previous results have indicated that these contaminants include various truncated, histidine-tagged forms of the target protein (data not shown). Accordingly, a second purification step was performed.

Purification of a histidine-tagged hydrolase expressed in *Pichia pastoris*


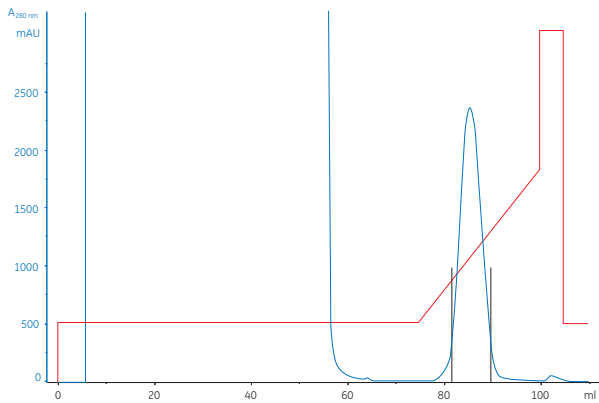


Fig 5. Black lines in chromatogram indicate the final pool. Arrow indicates the target protein band in the *Pichia* extract.

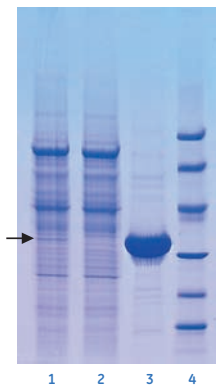

Lane 1: *Pichia* extract
Lane 2: Flow-through
Lane 3: Final pool
Lane 4: Low molecular weight marker

IMAC: 50 ml cell extract (corresponding to 20 g cells) applied to a 1 ml column with Ni Sepharose High Performance. 85 mM imidazole during equilibration, sample application and wash (see Optimization in Materials and Methods). Elution with a 25 ml linear imidazole gradient to 300 mM.

Results

- High-purity target protein was obtained from a low-expression sample.
- The Ni^{2+} concentration was low in the final pool of purified protein, 155 ppb (molar Ni^{2+} /protein ratio = 0.08).

Purification of 60 mg histidine-tagged protein on a 1 ml column

Comparison at high-load conditions between three different IMAC media


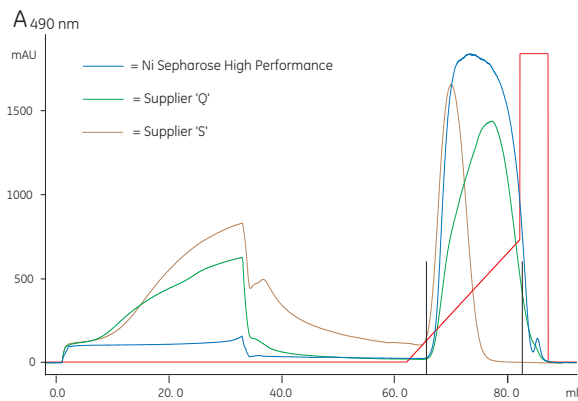


Fig 7. GFP-(His)₆ monitored at 490 nm. Black lines in chromatogram indicate the final pools.

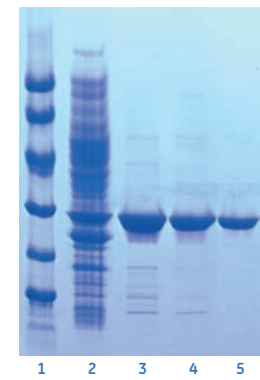

Lane 1: Low molecular weight marker
Lane 2: *E. coli* extract with GFP-(His)₆
Lane 3: Eluted from Ni sepharose High Performance
Lane 4: Eluted from Supplier 'Q'
Lane 5: Eluted from Supplier 'S'

IMAC: 32 ml extract applied to each 1 ml column. 5 mM imidazole during equilibration, sample application and wash. Elution with a 25 ml linear imidazole gradient to 250 mM. Pools of 16 ml were analysed.

Results

- Due to the high dynamic binding capacity, the yield of eluted protein for Ni Sepharose High Performance was almost twice that from the Supplier 'Q' medium, and over three times higher than that from the Supplier 'S' medium.
 - SDS-PAGE analysis of the final pools confirmed the superior binding capacity of Ni Sepharose High Performance.
- Note:* Improved purity can be achieved with a higher imidazole concentration than the 5 mM used in here binding buffer and sample.

Acknowledgements

We thank the owners of the clones used in this work:

Phadia, Uppsala, Sweden (MBP-(His)₆),

Dr. David Drew, Dept. of Biochemistry and Biophysics,
Stockholm University (GFP-(His)₆) and Dr. Mikael Widersten,
Protein engineering and redesign, Dept. of Biochemistry,
Uppsala University (*Saccharomyces hydrolase*).

www.gelifescience.com/hitrap

www.gelifesciences.com/protein-purification

GE Healthcare Bio-Sciences AB

Björkgatan 30

751 84 Uppsala

Sweden

GE, imagination at work, and GE monogram are trademarks of General Electric Company.

ÅKTAexplorer, Drop Design, ExcelGel, HiTrap, HisTrap, Sepharose and Superdex are trademarks of GE Healthcare companies.

All third party trademarks are the property of their respective owners.

Purification and preparation of fusion proteins and affinity peptides comprising at least two adjacent histidine residues may require a license under US pat 5,284,933 and US pat 5,310,663, including corresponding foreign patents (assignee: Hoffman La Roche, Inc).

© 2004-2007 General Electric Company - All rights reserved.
First published 2004

All goods and services are sold subject to the terms and conditions of sale of the company within GE Healthcare which supplies them. A copy of these terms and conditions is available on request. Contact your local GE Healthcare representative for the most current information.

GE Healthcare Europe GmbH
Munzinger Strasse 5, D-79111 Freiburg, Germany

GE Healthcare UK Ltd
Amersham Place, Little Chalfont, Buckinghamshire, HP7 9NA, UK

GE Healthcare Bio-Sciences Corp
800 Centennial Avenue, P.O. Box 1327, Piscataway,
NJ 08855-1327, USA

GE Healthcare Bio-Sciences KK
Sanken Bldg. 3-25-1, Hyakunincho, Shinjuku-ku,
Tokyo 169-0073, Japan

Asia Pacific T +85 65 62751830 F +85 65 62751829 • Australasia T +61 2 8820 8299 F +61 2 8820 8200 • Austria T 01 /57606 1613 F 01 /57606 1614 • Belgium T 0800 73 890 F 02 416 8206 • Canada T 1 800 463 5800 F 1 800 567 1008 • Central & East Europe T +43 1 972 720 F +43 1 972 722 750 • Denmark T +45 70 25 24 50 F +45 45 16 2424 • Eire T 1 800 709992 F +44 1494 542010 • Finland & Baltics T +358 9 512 3940 F +358 9 512 39439 • France T 01 69 35 67 00 F 01 69 41 98 77 Germany T 0800 9080 711 F 0800 9080 712 • Greater China T +852 2100 6300 F +852 2100 6338 • Italy T 02 26001 320 F 02 26001 399 • Japan T 81 3 5331 9336 F 81 3 5331 9370 • Korea T 82 2 6201 3700 F 82 2 6201 3803 • Latin America T +55 11 3933 7300 F +55 11 3933 7304 • Middle East & Africa T +30 210 96 00 687 F +30 210 96 00 693 • Netherlands T 0800-82 82 82 1 F 0800-82 82 82 4 • Norway T +47 815 65 777 F +47 815 65 666 • Portugal T 21 417 7035 F 21 417 3184 • Russia, CIS & NIS T +7 495 956 5177 F +7 495 956 5176 Spain T 902 11 72 65 F 935 94 49 65 • Sweden T 018 612 1900 F 018 612 1910 • Switzerland T 0848 8028 10 F 0848 8028 11 • UK T 0800 515 313 F 0800 616 927 • USA T +1 800 526 3593 F +1 877 295 8102



imagination at work