

SDS-PAGE Clean-Up Kit

Procedure A for sample volumes of 1 to 100 µL (containing 1–100 µg protein per sample)

Product protocol card

Introduction

Product codes

80-6484-70

Important

Read these instructions carefully before using the products.

Intended use

The products are intended for research use only, and shall not be used in any clinical or *in vitro* procedures for diagnostic purposes.

Description

Prior to starting procedure: Chill wash buffer (3) at -20°C for at least one hour. Add reductant to SDS-PAGE sample buffer (7) either 3.1 mg DTT or 5 µL 2-mercaptoethanol per 100 µL.

Step	Duration	Actions
1	15 min	Add 300 µL precipitant (1) to 1–100 µL sample (containing 1–100 µg protein per sample). Vortex. Incubate on ice. Add 300 µL co-precipitant (2) and mix.
2	5 min	Centrifuge (at least 12 000 × g). Tip: <i>Position tubes with cap hinge facing outward. Remove supernatant. Centrifuge briefly to bring remaining supernatant to bottom of tube. Remove remaining supernatant with micropipette.</i> Tip: <i>Reposition tubes as before. Proceed rapidly to avoid resuspension or dispersion of pellet. There should be no visible liquid remaining.</i>
3	5–10 s	Add 25 µL of water to the pellet. Vortex to disperse. Tip: <i>The pellet should disperse but not dissolve</i>
4	At least 30 min	Add 1 mL chilled wash buffer (3) and 5 µL of wash additive (4). Vortex 20–30 s every 10 min. Tip: <i>Make sure wash buffer is at -20°C before starting. The protein will not dissolve.</i>
5	5 min	Centrifuge (at least 12 000 × g).

6 No more than 5 min

Tip:

Position tubes with cap hinge facing outward. Remove and discard supernatant. Allow pellet to dry.

Tip:

Do not over-dry pellet.

7 5 min

Resuspend pellet in 5–40 µL buffer I (5). Vortex and incubate on ice.

8 5–10 min

Add 1 µL buffer II (6) for each 5 µL buffer I used in the previous step. Vortex and incubate on ice.

9 5–10 min

Add an equal volume (6–48 µL) SDS-PAGE sample buffer (7) with reductant. Vortex and incubate at room temperature

Tip:

If the solution turns yellow, add buffer I in 0.5 µL increments until the solution turns blue.

10 3 min

Heat at 95–100°C

Centrifuge briefly and gently tap tube to mix contents.

Sample is now ready to load on gel

cytiva.com

Cytiva and the Drop logo are trademarks of Global Life Sciences IP Holdco LLC or an affiliate.

All other third-party trademarks are the property of their respective owners.

© 2020 – 2021 Cytiva

All goods and services are sold subject to the terms and conditions of sale of the supplying company operating within the Cytiva business. A copy of those terms and conditions is available on request. Contact your local Cytiva representative for the most current information.

For local office contact information, visit cytiva.com/contact

29700228 AA V:4 05/2021

